



PRINCIPLE AND GENERAL DESCRIPTION OF THE MOUSE 4MID[®] KIT (Ref. 4VDX-18K8BB)

The Mouse 4MID[®] Kit (4VDX-18K3BB) is a sandwich ELISA assay that aims to detect and quantify proAKAP4 in murine sperm samples (in ejaculate, chilled or frozen semen in extender or in isolated spermatozoa). The Mouse 4MID[®] Kit in the BB format (4VDX-18K3BB) is composed of a 96-well microplate (12 x 8-wells strips), and all the reagents and buffers required to run the assay. The BB format is designed to run 2 samples per strip and then reused up to 12 times (12 strips). Therefore, the BB format enables running from 1 sample to 24 semen samples. The proAKAP4 is a sperm-specific protein that is a marker of sperm quality and male fertility in main mammals (Peddinti et al. 2008; Delehedde et al. 2018; Ruelle et al. 2019; Sergeant et al. 2019; Delehedde et al. 2019; Carracedo et al. 2020; Ruelle et al. 2020; Griffin et al. 2020; Bastan and Akcay, 2021; Dordas-Perpinyà et al. 2022; Boersma et al. 2022). Spermatozoa without proAKAP4 / AKAP4 are abnormal, immotile, and infertile (Miki et al. 2002; Fang et al. 2019; Delehedde et al. 2019). Oxidative stress decreases proAKAP4 concentrations in spermatozoa (Nixon et al. 2019; Delehedde et al. 2019; Sergeant et al. 2020). Each ELISA assay requires two major steps as described below in Part A and Part B of the Mouse 4MID[®] Kit - BB procedure. The proAKAP4 protein should be first released from spermatozoa using the Mouse Spermatozoa Lysis Buffer and according to the types and/or sperm concentrations of the samples (ejaculate, semen in extenders, or isolated spermatozoa) as described in Part A of the Mouse 4MID[®] - BB procedure. Then in Part B of the procedure, each prepared Mouse semen sample is loaded onto a 96-well microplate where a first antibody coated onto the bottom of the microplate will capture the proAKAP4 protein contained in the samples and will then be identified using a second antibody that is covalently coupled to horseradish peroxidase (sandwich method). A Substrate Solution is then added to each well and different color levels (blue) will appear proportionally to the concentrations of the proAKAP4 present in each semen sample. The color reaction is then stopped by adding the Stop Solution and the color intensity (yellow) is measured by spectrophotometry (reading at 450 nm). A positive control is included in each Mouse 4MID[®] Kit -BB. Ready-to-use Standard Solutions are provided to run a Linear Standard Curve allowing the calculation of proAKAP4 concentrations in Mouse semen samples. To compare animals or ejaculates or doses, we always recommend expressing proAKAP4 concentrations in ng per 10 million spermatozoa (ng/10M spz). Therefore, the concentrations in the number of spermatozoa in the samples should be known before running the assay or be performed independently of the Mouse 4MID[®] - BB assay.

Always read the following instructions carefully before use

TABLE OF CONTENTS

TABLE OF CONTENTS	1
I. REAGENTS AND MATERIALS INCLUDED	2
II. REAGENTS AND MATERIALS REQUIRED - NOT INCLUDED	2
III. STORAGE INFORMATION.....	2
IV. GENERAL INSTRUCTIONS OF USE	2
A. Preparation of the Semen Samples.....	2
a) Ejaculate or sample with sperm concentration above 50 million spermatozoa per mL Erreur ! Signet non défini.	
b) Frozen semen in extenders (chilled or frozen samples) or any sample with sperm concentration comprised between 10 to 50 million spermatozoa per mL.....	Erreur ! Signet non défini.
c) Isolated spermatozoa.....	Erreur ! Signet non défini.
B. Dosage of proAKAP4 Concentrations using the Mouse 4MID [®] Kit - BB.....	3
C. Calculation of Results – ProAKAP4 Concentration Determination	5
V. PRACTICAL ADVICE AND CAUTIONS	6



VI. REFERENCES	7
VII. RELATED PRODUCTS.....	8

I. REAGENTS AND MATERIALS INCLUDED

- R1 - Capture Antibody Coated Microplate of 96-wells (12 x 8-well strips)
- R2 - 1 Bottle of 10x Washing Buffer Solution (30 mL)
- R3 - 1 Bottle of 1x Dilution Buffer (30 mL)
- R4 - 5 Vials of Ready-to-Use Standard Solution (1.5 mL per vial)
- R5 - 1 Bottle of 1x Mouse Spermatozoa Lysis Buffer (30 mL)
- R6 - 6 Vials of Detection Antibody (110 µL)
- R7 - 1 Bottle of Substrate Solution (11 mL)
- R8 - 1 Bottle 1x Stop Solution (6 mL)
- R9 - 1 Vial of Positive Control (1.5 mL)
- R10 - 2 Adhesive Plate Sealers
- R11 - User Manual Instructions

II. REAGENTS AND MATERIALS REQUIRED - NOT INCLUDED

- 96-well microplate reader measuring absorbance at 450 nm
- Vortex
- Horizontal orbital microplate shaker (300 rpm)
- Polypropylene microtubes of 1.5 mL (for sample preparation)
- One glass bottle (for R2 dilution)
- One polypropylene tube of 15 mL (for R6 dilution)
- Multichannel pipette of 300 µL
- Pipette tips
- Pipettes of 20 µL, 200 µL and 1000 µL
- Plastic reagent reservoirs
- Scissors, pen, aluminum foil
- Ultrapure or double deionized water

III. STORAGE INFORMATION

- The Mouse 4MID® Kit - BB should be stored at 4°C upon receipt.
- All reagents must be protected from intense light.

IV. GENERAL INSTRUCTIONS OF USE

- Before use, bring all reagents except the R6 vial at **room temperature (RT)** for at least **30 minutes before running the assay**. The R6 Detection Antibody vial should always be kept at 4°C.
- Verify the absence of crystals in the R2 and the R5 bottles. In presence of crystals, gently agitate the solution until all crystals are completely dissolved. Then R2 and R5 solutions should be kept at ambient temperature.
- The R1 96-well Plate (12 x 8-well strips) is in a reusable aluminum foil pouch. The microplate frame and unused strips can be placed back in the reusable foil pouch for later use.

A. Preparation of the Semen Samples



ProAKAP4 should first be extracted from spermatozoa flagellum using a specific R5 Mouse Spermatozoa Lysis Buffer.

Do not forget before any pipetting of semen sample (ejaculate, chilled, or frozen semen in extenders or isolated spermatozoa) to resuspend cells by gently shaking the tube containing the semen as spermatozoa cells will pellet by sedimentation.

Please note that before lysing spermatozoa cells, the semen should always be brought to ambient temperature but **never kept on ice**.

a) Epididymis spermatozoa

1. After isolation of mouse epididymis spermatozoa, proceed to sperm numeration.
2. In a 1.5 mL conic tube, pellet the mouse spermatozoa (1 000 000 spermatozoa is required) by centrifugation at 2000 x g (5000 rpm) for 10 min at room temperature.
3. Remove the supernatant and add 190 µL of the R5 Mouse Spermatozoa Lysis Buffer.
4. Carefully vortex **continuously** each tube for 1 min at maximum speed (3000 rpm or above). All tubes should be vortexed before step 5.
Remark: Improper vortexing will lead to negative or false quantification value.
5. Add 200 µL of R3 Dilution Buffer to each sample.
6. Vortex 1 rapidly at maximum speed.
7. Keep the lysed samples at ambient temperature (17°C – 25°C) not more than 6 hours before use. Do not put it on ice.

b) Frozen spermatozoa

Unfroze the semen sample at 37°C for not more than 10 minutes, resuspend spermatozoa by gently mixing the semen and proceed as for epididymis spermatozoa.

B. Dosage of proAKAP4 Concentrations using the Mouse 4MID[®] Kit - BB

1. Open carefully the reusable aluminum foil pouch containing the R1 Microplate using scissors.
2. Add 100 µL of each Standard Solution (vial n°R4-1 to vial n° R4-5) on the first 8-well strip **from A1 well to E1 well** of the R1 Microplate (see table below). Please note that the R4-5 Standard Solution is the **Negative Control**. Always use a new tip for each different solution.

Standard vial n°	R4-1	R4-2	R4-3	R4-4	R4-5
ng/mL of proAKAP4	75	37.5	18.75	9.4	0
Well Position	A1	B1	C1	D1	E1

Always follow this plate scheme if using the calculation datasheet provided by 4BioDx[®] services.

3. Add **100 µL of the R9 Positive Control** in the F1 well of the first strip of the R1 Microplate.
4. Then add **150 µL of one or two semen samples as prepared in Part A** in the G1 and H1.

Remark 1: Vortex rapidly each semen sample before loading

Remark 2: Always use a new tip for each different semen sample.

5. Cover the R1 Microplate with one R10 Adhesive Plate Sealer and incubate the R1 Microplate **for 2 hours** at room temperature on a horizontal shaker with gentle shaking (200 rpm).



6. During the incubation time:

- a. Prepare the Detection Antibody: Dilute the content of one R6 Detection Antibody tube in 990 μ L of R3 Dilution Buffer in a 1.5 mL tube. Close the tube and mix the Detection Antibody Solution by reversing the tube several times. Keep this solution at ambient temperature until Step 9.

Remark 1: Before dilution, please carefully hit the bottom of the R6 tube vertically or rapidly centrifuge to pull down any liquid in the cap of the R6 tube. One tube can be used for two strips / two assays.

Remark 2: After dilution in R3, the Detection Antibody Solution should be used within one month and kept at 4°C.

- b. Prepare the Washing Solution: Add 30 mL of the R2 into 270 mL of ultrapure water (or double deionized water) in a clean glass bottle. Gently agitate the solution, avoid foaming and keep it at ambient temperature.

Remark: After use, the remaining Washing Solution should be kept at 4°C between each use.

7. At the end of incubation time, remove the R10 Adhesive Plate Sealer and eliminate the sample solutions by reversion of the R1 Microplate (or aspirate when using an automatic microplate washer).

8. Transfer some Washing Buffer to a plastic reagent reservoir. Wash each well by adding **300 μ L of R2 Washing Buffer 1x Solution** using the multichannel micropipette. Then discard the Washing Solution by reversion of the R1 Microplate. Repeat two times more. Please tap down gently the R1 Microplate on an absorbent dry paper to remove residual R2 Washing Buffer droplets between each washing step and before loading the R6 Detection Antibody.

Remark: Results of the assay will markedly be influenced by the proper performance of the washing steps.

9. Transfer all the Detection Antibody Solution to a plastic reagent reservoir. Add **100 μ L of Detection Antibody Solution** prepared in step 6 to each 96-well Microplate (R1) well using the multichannel micropipette.

10. Cover the R1 Microplate with a new R10 Adhesive Plate Sealer and incubate the R1 Microplate **for one hour** at room temperature with gentle shaking (300 rpm).

11. Remove the R10 Adhesive Plate Sealer and eliminate the Antibody Solution by reversion of the R1 Microplate (or aspiration when using an automatic microplate washer).

12. Transfer some Washing Buffer to a plastic reagent reservoir. Wash each well by adding **300 μ L of R2 Washing Buffer 1x Solution** using the multichannel micropipette. Then discard the Washing Solution. Repeat two times more. Please tap down gently the R1 Microplate on an absorbent dry paper to remove the residual Washing Buffer droplet between each washing step and before adding the R7 Substrate.

Remark: Results of the assay will markedly be influenced by the proper performance of the washing steps.

13. Transfer the Substrate Solution to a plastic reagent reservoir. Add **100 μ L of R7 Substrate Solution** to each well using the multichannel pipette.

Remark 1: Keep carefully away from light with an aluminum sheet.



Remark 2: Please note that R7 Substrate Solution and R8 Stop Solution (step 15) should be added to wells in the same order and with the same time interval.

- 14.** Protect from light and incubate the R1 Microplate under gentle shaking (300 rpm) for **10 minutes** at room temperature.

Remark: The incubation must not exceed 30 minutes.

- 15.** Transfer the Stop Solution to a plastic reagent reservoir. Add **50 µL of R8 Stop Solution** to each well using the multichannel pipette and place the R1 Microplate for 1 minute on an orbital shaker at 300 rpm to mix well before reading the R1 Microplate.

- 16.** Determine the samples' optical density using a microplate reader set to 450 nm.

*Remark: Always perform the measure **immediately** after adding the Stop Solution.*

C. Calculation of Results – ProAKAP4 Concentration Determination


A Standard Curve must be performed for each analysis to calculate proAKAP4 concentrations in Mouse semen. You can either do the Standard Curve a) using the Calculation Data Sheet provided for each sperm preparation protocol (part A) or b) do it manually yourself using software such as Excel.

Remark: always take into account the dilutions you have made when considering sperm concentrations of the semen samples.

a) Using the Calculation Data Sheets provide by 4BioDx® services (upon request at contact@4biodx.com). Each Calculation Data Sheet has been designed to automatically draw the Standard Curve from your optical density values (OD) and calculate the proAKAP4 concentrations first in ng /mL, then in ng/10 million spermatozoa. First select the calculation data sheet according to the sample preparation procedure (a, b, or c) you have used in Part A. You will then need to fill the blue cases with OD obtained from the plate reader and follow the indicated steps on the sheet to draw the curve and then the sheet will automatically calculate the concentrations of proAKAP4 in ng/mL. By adding the concentration of spermatozoa in million per mL of the semen samples, or the number of spermatozoa in the case of isolated spermatozoa (blue column of step 5), you will obtain the concentrations of proAKAP4 in ng/10 million spermatozoa. You may then copy and paste the values, tables, and graphs into your own report files. Please note that hitting the step 6 button will generate a new sheet with the summary of the results.

b) Manually, you will need to first subtract the optical density obtained from value E1 to the optical density of each standard (A1 to E1) and each sample (G1 and H1) that you obtained from the plate reader. Then create a Standard Curve by reporting the data (optical density values) on an Excel Spreadsheet. Express the optical densities (OD) in abscissa in the function of the proAKAP4 quantities in ng/mL in ordinate and then generate a two-degree polynomial regression equation (ProAKAP4 [conc] = a x (OD)² + b x (OD)) using the data analysis tools of an Excel Spreadsheet.

ProAKAP4 in (ng/mL)	Optical density (OD) values
75	OD A1
37.5	OD B1
18.75	OD C1
9.4	OD D1
0	OD E1

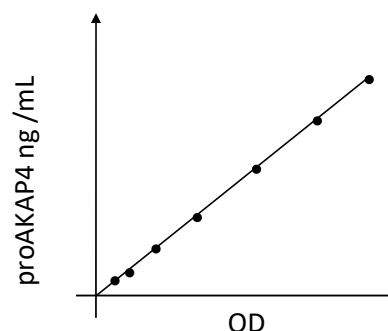



Then to **express the proAKAP4 concentrations in ng/mL**, calculate the concentrations using this two-degree polynomial regression equation.

To compare animals, ejaculates, or any semen sample in extenders, we recommend expressing proAKAP4 concentrations **in ng per 10 million spermatozoa (ng/10 M spz)**.

To express results in ng of proAKAP4 per 10 million spermatozoa, the formula will be:
 Conc. proAKAP4 ng/10M = (conc. proAKAP4 in ng/mL / conc. spz M/mL) x 10 x dilution factor* x (2/3)
 (*Use the dilution factor of 13.3 and 10 for the sample preparation procedure a) or b), respectively).

In the case of isolated spermatozoa: you should first calculate the concentration of spermatozoa in Million per mL (M/mL). To do so, multiply by 2.5 the spermatozoa number contained in the pellet, normally containing 1 000 000 spermatozoa at step 2a of the Semen sample preparation part.

Express then the results in ng of proAKAP4 per 10 million spermatozoa (ng/10M spz) with the following formula:
 Conc. proAKAP4 in ng/10M of spz = (conc. proAKAP4 in ng/mL / conc. spz M/mL) x 10 x 6.66

Example of threshold values of proAKAP4 concentrations in rodent semen (all breeds):

ProAKAP4 concentrations	Semen Quality	Long Lasting Motility
Less than 15 ng/10M spz	Poor	+/-
Between 15 and 40 ng/10M spz	Good	+
Between 40 and 60 ng/10M spz	Very Good	++
Over 60 ng/10M of spz	Excellent	+++

V. PRACTICAL ADVICE AND CAUTIONS

- For research use only purposes and not for use for diagnostic purposes.
- Always follow good laboratory practices.
- Always follow the plating scheme indicated in part B when using the Calculation Datasheet provided by 4BioDx services.
- Always use a new tip for each different solution and for each semen sample.
- The use of a multichannel pipette is mandatory to ensure the timely delivery of liquids.
- Use the supplied reagents as an integral unit prior to the expiration date.
- After dilution in R3, the Detection Antibody Solution should be used within one month and always kept at 4°C.
- To avoid distortions due to differences in incubation times, R7 Substrate Solution and R8 Stop Solution should be added to wells in the same order and with the same time interval.
- The R7 Substrate Solution can be irritating to the skin.
- The R8 Stop Solution can be harmful in case of ingestion and could lead to irritation when in contact with the skin.
- Do not expose the Substrate Solution to light or oxidative substances.
- Use only reagents from the same Mouse 4MID® Kit - BB.
- Any variation in ambient temperature, pipetting, washing method, or incubation time can cause variation in optical density results.
- Observe all federal, state, and local regulations in terms of waste disposal.



VI. REFERENCES

- Boersma A, Primus J, Wagner B, Broukal V, Andersen L, Pachner B, Dahlhoff M, Rüllicke T and Auer KE (2022)** Influence of sperm cryopreservation on sperm motility and proAKAP4 concentration in mice. *Reprod Med Biol.* Vol. 21:e12480. doi: 10.1002/rmb2.12480
- Dordas-Perpinyà M, Sergeant N, Ruelle I, Bruyas JF, Charreaux C, Michaud S, Carracedo S, Catalán J, Miró J, Delehedde M and Briand-Amirat L (2022)** ProAKAP4 semen concentrations as a valuable marker protein of post-thawed semen quality and bull fertility: a retrospective study. *Veterinary Sciences.* Vol. 9 :224.
- Kowalczyk A, Gałeska E and Bubel A (2022)** The concentration of proAKAP4 and other indicators of cryopotential of spermatozoa cryopreserved in extender with holothuroidea extract addition. *Animals.* Vol. 12:521.
- Riesco MF, Neila-Montero M, Palacín-Martínez C, Montes-Garrido R, Álvarez R; Boixo JC; De Paz P and Anel L and Anel-López L (2022)** Establishment of innovative biomarkers to optimize cooling and cryopreservation protocols in ram sperm. *Reproduction in Domestic Animals.* Vol. 57(1):101.
- Dordas-Perpinya M, Yanez I, Catalan J, Eddarkaoui S, Delehedde M, Sergeant N, Bruyas JF, Briand-Amirat L, and Miró J (2022)** ProAKAP4 concentrations in stallion and donkey: Comparison with kinetic parameters, motility and concentration. *Reproduction in Domestic Animals.* Vol. 57(1):61.
- Malo C, Carracedo S, Delehedde M, Sergeant N and Skidmore L (2021)** Identification of proAKAP4 in Dromedary sperm and their correlation with monthly semen parameters. *Reproduction & Fertility.* Vol.2(4):268-279.
- Bastan I and Akcay E (2021)** Quality assessment of frozen bull semen with the precursor A-kinase anchor protein 4 biomarker. *Andrologia* e14164.
- Blommaert D, Sergeant N, Delehedde M, Donnay I, Lejeune JP, Franck T and Serteyn D (2021)** First results about ProAKAP4 concentration in stallion semen after cryopreservation in two different freezing media. *Cryobiology.* Vol. 102:133-135
- Carracedo S, Loyens A, Eddarkaoui S, Serteyn D, Malo C, Skidmore L, Briand-Amirat L, Barbotin AL, Maurage CA, Delehedde M and Sergeant N (2020)** The sperm specific proAKAP4 polypeptide exhibited conserved functions, localizations and metabolism among mammals. *Animal Reproduction Science.* Vol. 220 :106448
- Riesco M, Anel-Lopez L, Neila-Montero M, Palacin-Martinez C, Montes-Garrido R, Alvarez M, de Paz P, Anel L (2020)** ProAKAP4 as Novel Molecular Marker of Sperm Quality in Ram: An Integrative Study in Fresh, Cooled and Cryopreserved Sperm. *Biomolecules.* 10(7):1046.
- Wu S, Mipam T, Xu C, Zhao W, Shah MA, Yi C, Luo H, Cai X, Zhong J (2020)** Testis transcriptome profiling identified genes involved in spermatogenic arrest of cattleyak. *PLoS One.* 15(2):e0229503.
- Ruelle I, Sergeant N, Bencharif D, Charreaux F, Thorin C, Michaud S, Dordas-Perpinyà M, Jouy N, Audry S, Maurage CA, Delehedde M and Briand-Amirat L (2020)** ProAKAP4 concentrations in semen as a predictive tool of bull fertility: A preliminary study. *Reproduction, Fertility and Development.* 32(1):145.
- Sergeant N, Carracedo S, Blommaert D, Aubry S, Jouy N, Lejeune JP, Franck T, Maurage CA, Serteyn D, Buée L and Delehedde M (2020)** Proteolysis of proAKAP4 in semen as a regulatory sensor of sperm quality and functionality. *Andrology.* Vol. 8 (Suppl. 1):44-45.
- Nixon B, Bernstein I, Cafe SL, Delehedde M, Sergeant S, Eamens AL, Lord T, Dun MD, De Iuliis GN and Bromfield EG (2019b)** A Kinase Anchor Protein 4 is vulnerable to oxidative adduction in male germ cells. *Frontiers in Cell and Developmental Biology* – Vol. 7:319.
- Fang X, Huang LL, Xu J, Ma CQ, Chen ZH, Zhang Z, Liao CH, Zheng SX, Huang P, Xu WM, Li N, Sun L (2019)** Proteomics and single-cell RNA analysis of Akap4-knockout mice model confirm indispensable role of Akap4 in spermatogenesis. *Developmental Biology.* 454(2):118-127.
- Sergeant N, Briand-Amirat L, Bencharif D and Delehedde M (2019)** The sperm specific protein proAKAP4 as an innovative marker to evaluate sperm quality and fertility. *Journal of Dairy & Veterinary Sciences.* Vol. 11:01-19.



Singh R, Junghare V, Hazra S, Singh U, Sengar GS, Raja TV, Kumar S, Tyagi S, Das A.K., Kumar A, Koringa P, Jakhesara S, Joshi CJ and Deb R (2019) Database on spermatozoa transcriptogram of catagorised frieswal crossbred (holstein friesian X sahiwal) bulls. *Theriogenology*. Vol. 129:130-145.

Ruelle I, Charreaux F, Bencharif D, Thorin C, Michaud S, Schmitt E, Sergeant N, Briand-Amirad L and Delehedde M (2019) Assessment of the sperm specific protein proAKAP4 as a marker to evaluate sperm quality and fertility in Holstein bulls. *Rev. Bras. Reprod. Anim.* Vol.43(2):472.

Delehedde M, Carracedo S, Selleslagh M, Eddarkaoui S, Amirat-Briand L and Sergeant N (2019) ProAKAP4 polypeptide as a biomarker of sperm functionality and male fertility disorders. *Int J Gynecol and Reprod Sci.* Vol. 2(1):13-19.

Delehedde M, Demouveaux B, Remy G, Selleslagh M, Dewulf Q, Desseyn JL, Moreau M, Gosset P, Muriel Pichavant M and Sergeant N (2019) ProAKAP4 concentrations as an indicator of good spermatogenesis and sperm quality under oxidative stress conditions. *Andrology*. Vol. 7(1):86.

Fu Q, Pan L, Huang D, Wang Z, Hou Z and Zhang M (2019) Proteomic profiles of buffalo spermatozoa and seminal plasma. *Theriogenology*. Vol. 134:74-82.

Delehedde M, Blommaert D, Jouy N, Scabello J, Miersman H, Franck T, Sertejn D, Mitchell V, and Sergeant S (2018) Concentration of proAKAP4 as a pertinent read-out of sperm quality in mammals. *Animal Reproduction Science*. Vol. 194:24.

Miki K, Goulding EH, Fulcher KD, Eddy E, Willis WD and Brown PR (2002) Targeted disruption of the AKAP4 gene causes defects in sperm flagellum and motility. *Developmental Biology*. Vol. 248:331-342.

VII. RELATED PRODUCTS

Reference	Designation	Specificity
4VDX-18K2	Pig 4MID [®] Kit	Pig proAKAP4
4VDX-5 / 4VDX-18K3BB	Horse 4MID [®] Kit	Horse / Donkey proAKAP4
4VDX-18K4	Bull 4MID [®] Kit	Bull proAKAP4
4VDX-18K5 / 4VDX-18K5BB	Dog 4MID [®] Kit	Dog proAKAP4
4VDX-18K6	Rabbit 4MID [®] Kit	Rabbit proAKAP4
4VDX-18K7	Ram 4MID [®] Kit	Ram proAKAP4
4BDX-18K8 / 4BDX-18K8BB	Mouse 4MID [®] Kit	Mouse / Rat proAKAP4
4VDX-18K9	Goat 4MID [®] Kit	Caprine proAKAP4
4VDX-19K10	Cat 4MID [®] Kit	Cat proAKAP4 (feline)
4VDX-19K11	Camel 4MID [®] Kit	Camelids proAKAP4

FOR RESEARCH USE ONLY PRODUCT / NOT FOR USE IN DIAGNOSTIC PROCEDURES