FOOD ALLERGEN SCREENING TEST KIT

FAST KIT ELISA VER. III





THIS PRODUCT CAN MEASURE "EGG", "PEANUT",
"WHEAT", "BUCK WHEAT", "SOYBEAN", "MILK" AND
"SESAME" PROTEINS IN FOOD SUCH AS RAW
MATERIALS AND PROCESSED FOODS WIDELY.

FEATURES

- High precision quantity measurement is possible.
- Kit has been improved by removal of 2-mercaptoethanol from "FASTKIT ELISA Ver. II.
- Equal examination operation as "FASTKIT ELISA Ver. II"



FAST KIT ELISA VER.III

FOOD ALLERGY MATERIAL SCREENING EXAMINATION KIT

An example of measurement from food

1g of crushed food sample

Add 19ml of Extraction medium and concuss it more than 12 hours

After centrifugation, remove inessential materials by filtration

Dilute it to 20 times with dilution buffer

ELISA operation

Measure absorbance with plate leader

Description	Cat. No.
FASTKIT ELISA Ver. III EGG	NPH-999100430EX
FASTKIT ELISA Ver. III WHEAT	NPH-999100431EX
FASTKIT ELISA Ver. III PEANUT	NPH-999100432EX
FASTKIT ELISA Ver. III BUCKWHEAT	NPH-999100433EX
FASTKIT ELISA Ver. III MILK	NPH-999100434EX
FASTKIT ELISA Ver. III SOYBEAN	NPH-999100435EX
FASTKIT ELISA Ver. III SESAME	NPH-999100436EX

Quantity: 1 kit (96 wells (8 wells ×12 strips))

Storage: 2-8℃


This product is based on the guidelines shown by CAA Notification No. 286 from the Deputy Secretary General of the Consumer Affairs Agency dated September 10, 2010, "Test methods for food products containing allergic substances".

FASTKIT SLIM SERIES



Anybody can perform the test easily!

Applicable for smear tests!

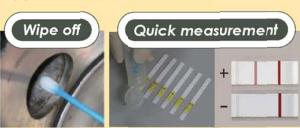

Detect allergy substances in 15 minutes!

Product Features

- Simple one-step operation that only requires a drop of sample solution
- Judgment: appearance of reddish violet line
- Simple and rapid test is best for routine management at the production sites.

Description	Cat. No.
FASTKIT SLIM EGG	NPH-NFS001
FASTKIT SLIM MILK	NPH-NFS002
FASTKIT SLIM WHEAT	NPH-NFS003
FASTKIT SLIM BUCKWHEAT	NPH-NFS004
FASTKIT SLIM PEANUT	NPH-NFS005
FASTKIT SLIM SOYBEAN	NPH-NFS006

Quantity: 20 tests Storage: 2-8°C

6

COSMO BIO CO., LTD.

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME, KOTO-KU. TOKYO 135-0016, JAPAN

TEL: (81)3-5632-9617
FAX: (81)3-5632-9618
e-mail: export@cosmobio.co.jp
URL: www.cosmobio.com



10122