*For research use only. Not for clinical diagnosis.*

**Catalog No. CK021**

**Basic Medium for Porcine Oocyte Maturation (POM)**

**BACKGROUND**

This is a modification medium which added polyvinyl alcohol to the original medium for porcine oocyte maturation (POM) (Yoshioka et al. <sup>1)</sup>) and is a complete synthetic basic medium for oocyte maturation. This medium effectively supports porcine oocyte maturation under a low oxygen condition (5%O<sub>2</sub>/5%CO<sub>2</sub>/90%N<sub>2</sub>, 39°C).

- Usage** Please add 10 IU/ml equine chorionic gonadotropin (eCG) + 10 IU/ml human chorionic gonadotropin (hCG) to the POM medium during porcine oocyte maturation culture. There is no need to add porcine follicular fluid.
- POM Composition** Mineral salts, Sugars, Sodium bicarbonate, Amino acids, Hypotaurine, polyvinyl alcohol , etc.
- Storage** Stored at 4°C in a dark room. Do not freeze.
- Expiration date** The expiration date is shown on the label of the medium bottle (4 months after the medium preparation)
- Package** 25 ml × 3 bottles
- Note** POM medium is for research use only. Do not use in animals, humans, or diagnostic procedures.
- References** 1) Yoshioka. K, Suzuki. C, Onishi. A: J. Reprod. Dev., 54: 208-213, 2008.

*For research use only. Not for clinical diagnosis.*

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COSMO BIO Co., LTD.