For research use only. Not for clinical diagnosis.

Catalog No. CK022

High Performance Basic Medium for Porcine Oocyte Maturation (HP-POM)

BACKGROUND

HP-POM is a high-performance medium that added TGF- α to POM (#CK021) and is a serum-free basic medium for porcine oocyte maturation. Although the rate of oocyte maturation in this medium is almost at equal as POM, the increased rate of embryonic development after in vitro fertilization (the blastocyst formation) is reported ¹⁾.

This medium is sterilized and effectively supports porcine oocyte maturation under a low oxygen condition (5%O₂/5%CO₂/90%N₂, 39°C).

Usage	Please add 10 IU/ml equine chorionic gonadotropin (eCG) + 10 IU/ml human chorionic gonadotropin (hCG) to HP-POM medium during porcine oocyte maturation culture. There is no need to add porcine follicular fluid.
POM Composition	Mineral salts, Sugars, Sodium bicarbonate, Amino acids, Hypotaurine, polyvinyl alcohol, TGF- α etc.
Storage	Stored at 4°C in a dark room. Do not freeze.
Package	25 ml × 3 bottles
Expiration date	The expiration date is shown on the label of the medium bottle (4 months after the medium preparation)
Note	HP-POM medium is for research use only. Do not use in animals, humans, or diagnostic procedures.
References	1) Yoshioka. K, Suzuki. C, Onishi. A: J. Reprod. Dev., 54: 208-213, 2008.

For research use only. Not for clinical diagnosis.

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COSMO BIO Co., LTD.