for **IMMUNOSTAINING**

A-PAP PEN

ADVANCED PAP PEN

PAP PEN minimize the consumtion of liquid for immunostaining by using this PAP PEN as a "fence" on slide glass.

It is very easy to use. You only write a "circle" around tissue by this pen on slide glass. Then, a fence are created on the slide glass to prevent overflow of liquid for immunostaining.

Usage

Before fist use:

- 1. Before first use, hold the pen tip upside and press the tip down for air vent.
- 2. Hold the pen upright to press the tip down only once until the tip becomes saturated.
- 3. Do 1 again to reject ink on the tip holder, and wipe overflowing ink around the tip.
- 4. Remove the ink that has just come out first time.
- * Be sure to close the cap after use.

Two pen tip sizes, 4mm(Regular) and 2mm (Mini).

REACH compliant

T

Precautions for use

- · When using for the first time, follow the instructions on the pen body to bleed air before using.
- · Please perform the dyeing work after the ink has dried well. If too much ink comes out, wipe it off.
- It will not function if the surfactant remains on the glass after immersion in PBS.

Wipe it off or immerse the glass in a buffer solution that does not contain a surfactant and let the surfactant flow.



