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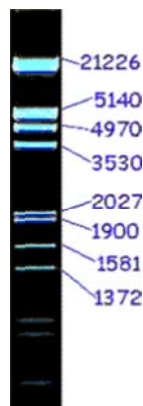
## Bio-Imager Pen

Catalog No. DAI-BIP

The Bio-Imager Pen, which contains a nontoxic, phosphorescent writing material is designed to record data and position information on membranes prior to exposure to X-ray film, allowing permanent records. The pen has widespread use in molecular biology reducing waste, saving time and avoids the use of isotopic materials. While the writing material casts a greenish white glow in the dark, it is recharged by light.

### Principle

Letters or symbols written using the Bio-Imager Pen absorb light when in light conditions and release the energy absorbed as light when in dark conditions. Film is exposed to this light and is developed as an image.

### Application

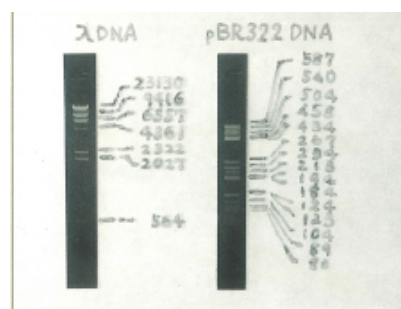
Use the pen to record letters or symbols directly on the DNA ladder or on cling wrap covering the DNA markers.

### Directions to Use

To use the Bio-Imager pen, first shake the pen horizontally, and then press the tip straight down several times until the fluid starts soaking the tip. The pen is then ready for use. Allow the writing to dry for approximately 1-2 minutes at room temperature, and then expose to the X-ray film. Optimal images are produced in approximately 10-12 hours.

### Precaution

When not in use keep the pen tightly closed, and store in a horizontal position.

COSMO BIO Co., LTD.  
Inspiration for Life Science

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME, KOTO-KU, TOKYO 135-0016 JAPAN

[International/Outside Japan]

TEL: (81)3-5632-9617 / FAX: (81)3-5632-9618

e-mail: export@cosmobio.co.jp / URL: www.cosmobio.com

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TEL: 03-5632-9610 / FAX: 03-5632-9619

e-mail: support@cosmobio.co.jp / URL: www.cosmobio.co.jp