[Cell Freezing Medium] (Chemically defined)
[Cell Freezing Medium]

COS banker
COS banker II

Catalog No. KOJ-COS-CFM01

Catalog No. KOJ-COS-CFM02

COS banker is a one-step culture cell storage solution. Suspend culture cells in COS banker and store at -80°C for long term storage (it is not necessary to store in liquid nitrogen). This is a serum-free solution with no animal products added.

Features

- Ready-to-use
No dilutions required
Store in -80°C freezer
Rapid and long term storage at -80°C
No serum
No animal products




Instructions

- 1. Collect culture cells (during logarithmic growth stage) and place 5x10^5 - 1x10^7 cells in ... ge for 5 minutes at 1,200 rpm.
2. Add 1 ml of COS banker to cell pellet and vigorously mix, then freeze and store at -80°C.
3. If liquid Nitrogen storage is desired, freeze cells at -80°C first and then transfer to liquid Nitrogen.
4. To thaw, defrost in 37°C water bath quickly.

Note: It is ideal to use a logarithmic growth stage culture in order to increase the survival rate when thawed.

Contents and Usage

Table with 3 columns: Product Name, COS banker, COS banker II. Rows include Product Code, Size, Specification, Storage, Expiry Date, and Application Tested Cell.

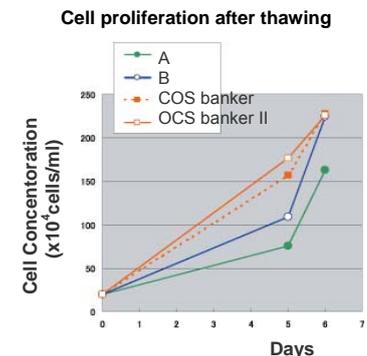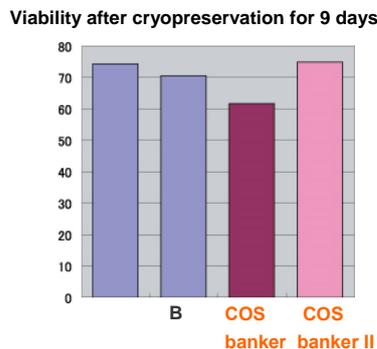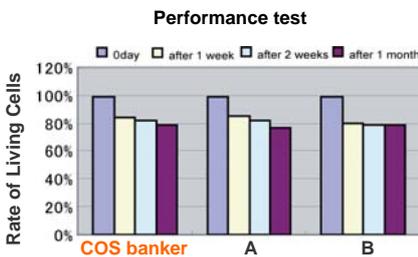Quality Controls

Endotoxin-free tested / Bacteria free tested

Warnings

- Research use only
Not for diagnostic and/or clinical use

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[Cell Freezing Medium] (Chemical defined)

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COSMO BIO CO., LTD.