**For research use only. Not for clinical diagnosis.**

**Catalog No.PMC-SWN-IMS32C**

# Mouse Schwann Cell Line (IMS32)

## Cell specification

Schwann cells are the principal glia cells of the peripheral nervous system (PNS) and are involved in several aspects of PNS biology. In addition to providing trophic support for neuronal development and maintenance through the production of cytokines and growth factors, Schwann cells wrap around axons of motor and sensory neurons to create an insulating myelin sheath that aids impulse conductivity. In recent years, Schwann cells have increasingly known for their support to nerve regeneration while their dysfunction may contribute to Neurodegenerative diseases such as ALS.

IMS32s are spontaneously immortalized cell line established from long term culture of mouse dorsal root ganglia and peripheral nerves. IMS32 cells retain many characteristics of mature Schwann cells and are readily propagated in the supplied media. IMS32 cells can be used to investigate many aspects of neuronal biology including nerve regeneration, neuroprotective mechanisms, and for the development of novel therapeutic approaches to neurological disorders.

## Components

Product Name	Quantity	Amount	Storage Conditions	Stability
Mouse Schwann Cell Line (IMS32), Cryopreserved	5 x 10 <sup>5</sup> cells/vial	1 vial	Liquid nitrogen storage	6months

- ※ Store frozen cells in liquid nitrogen if not ready for use after receipt.
- ※ Based on the license agreement of Tokyo Metropolitan Institute of Medical Science, IMS32 cell are prohibited to provide (distribution, lending, transfer, licensing, etc.) to a third party.

## General Information

<b>Organism</b>	ICR mouse (adult)
<b>Tissue</b>	Dorsal Root Ganglion and Peripheral Nervous System
<b>Cultural Properties</b>	Adherent
<b>Biosafety Level</b>	1
<b>Morphology</b>	Spindle-shaped
<b>Cell Type</b>	Spontaneous Immortalization
<b>Quality Check</b>	Mycoplasma (-), Immunohistochemical reactivity of cell markers

## Optimized Culture Medium (selling separately)

Catalog No.	Product Name	Size	Quantity	Storage Conditions	Expiration date
SWNMM	Schwann Cell (IMS32) Medium	500 mL	1	-20°C Freezer	- Written on the bottle
				4°C	3month after thawing

## Protocols

- ※ This product is available for passage.

### [Thawing Protocol]-for 100mm plastic culture dish

Preparation

- 100mm dish for culturing cells
- Medium for the mouse Schwann cell line (IMS32)

1. Thaw vials of frozen cells by warming them in 37°C warm water for 2 minutes.
2. Transfer the thawed cell solution to a 50 mL centrifuge tube containing 10 mL of medium previously returned to room temperature, mix, and collect the cell solution by 1 mL separating the culture medium in the centrifuge tube and co-washing the vials.
3. Transfer cell suspension to 100mm dish (seeding density: approximately  $0.9 \times 10^4$  cells/cm<sup>2</sup>). Incubate the plate at 37°C under 5% CO<sup>2</sup>

Note: Centrifugation of cells after thawing are not recommended since these actions are more harmful to the cells than the effect of residual cell freezing medium in the culture and cell attachment.

4. Replace the medium with medium heated to 37°C once the day after seeding, and then 2 to 3 times a week.

\*After 3~5 days, IMS32 cells reach approximately 80~100% confluent.

### **[Cell passage Protocol]**

Preparation

- 70-90% confluent cells
- Medium for the mouse Schwann cell line (IMS32)

1. Remove 70-90% confluent cells from the CO<sub>2</sub> incubator.
2. Aspirate the supernatant, add 10-12mL/100mm medium.
3. Detouch the cells by pipetting.  
※ After pipetting, check under a microscope to see if the cells are detached. If pipetting is inadequate, a significant number of cells cannot be detached, and growth may be affected.
4. Collect the cell suspension in a 50mL centrifuge tube.
5. Mix the cell suspension and medium in a ratio of 1:3 to 1:4 and seed it in a new culture vessel.
6. Incubate in a 37°C incubator in the presence of 5% CO<sub>2</sub>. Passage should be performed in the same manner with cells 70-90% confluent (see Figure 1, C).


### **[Freezing Cells Protocol]**

1. Cell passage protocols 1 to 4 are performed.
2. Centrifuge at 4°C for 5 minutes at 200 ×g, remove the supernatant, add COS banker(KOJ product cord COS-CFM01) to  $1 \times 10^6$  cells/mL and suspend.
3. Dispense 1 mL/tube into cryopreservation tubes, and cryopreserve at -80°C by using cell freezing containers (BM-instrument-product-code BCS-136 or equivalent products).
4. Store at -80°C for 1 day or more. Transfer to liquid nitrogen.


## Technical information


A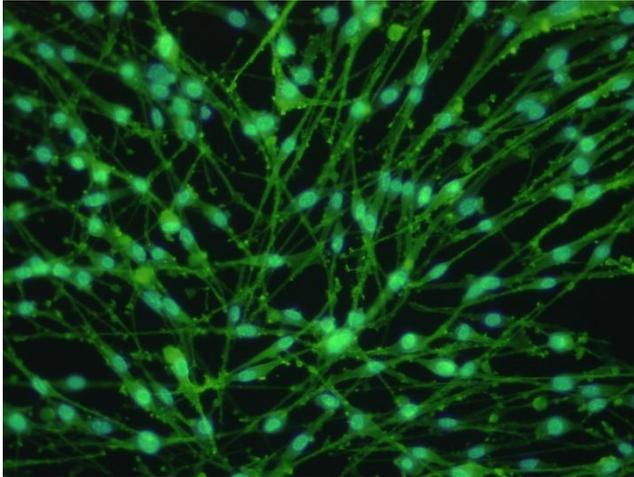

B

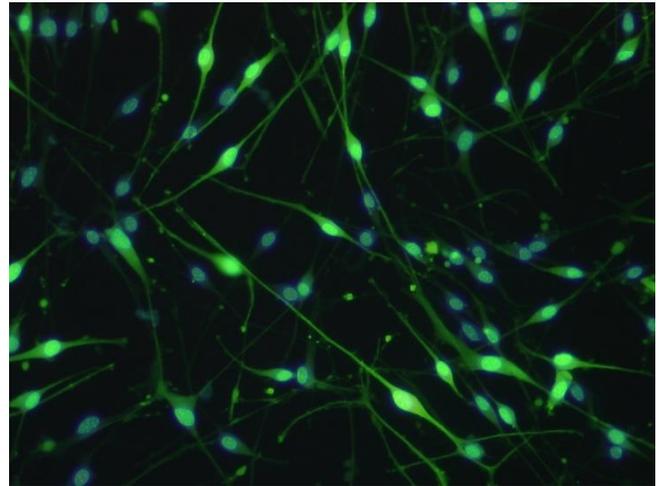C
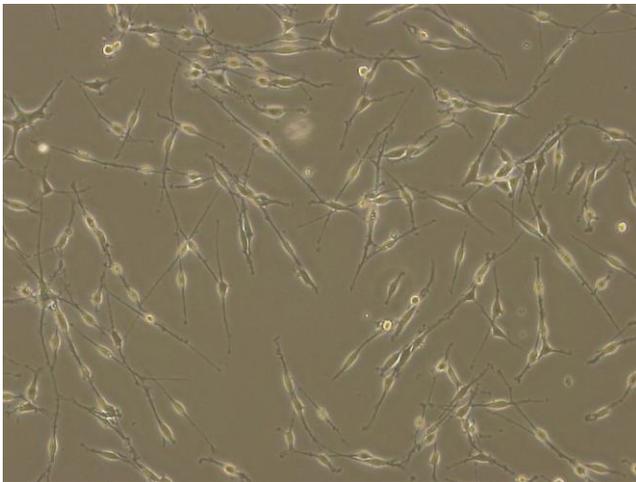
### [Immunofluorescence of IMS32 cells]

A: Anti-p75NTR (green), nuclear staining (blue)

B: Anti-s100 (green), nuclear staining (blue)

C: phase-contrast microscopic image

Magnification :  $\times 100$ (A,B,C)

## References

- 1) Watabe K, Fukuda T, Tanaka J, et al. Spontaneously immortalized adult mouse Schwann cells secrete autocrine and paracrine growth-promoting activities. *J Neurosci Res.* 1995;41:279-290.
- 2) Watabe K, Sakamoto T, Kawazoe Y, Michikawa M, Miyamoto K, Yamamura T, Saya H, Araki N. Tissue culture methods to study neurological disorders: Establishment of immortalized Schwann cells from murine disease models. *Neuropathology* 2003;23:64-74.

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COSMO BIO Co., LTD.