

Ab-Carrier™

Transfection with Anti-nuclear pore complex (NPC) antibody inhibits HeLa cell proliferation

Experimental protocol

Seeding HeLa cells in 96-well plate

(5×10^4 cells/mL; medium volume MEM (+ 10% FBS) 0.1 mL/well)

↓ Incubate at 37° C in the presence of 5% CO₂ for 24 hours

Ab-Carrier™ 0.5 μL was added to 10 μL of anti-NPC antibody (0.1 mg/mL) and mixed well.

↓ room temperature for 20 minutes

Add 10.5 μL/well of reaction solution to HeLa cells after 24-hour culture.

↓ 37° C in the presence of 5% CO₂ for 4 hours

Remove the medium and wash with PBS 0.1 mL/well × 2 times.

↓

Add 0.25% Trypsin-EDTA 20 μL/well

↓ 37° C in the presence of 5% CO₂ for 2 minutes

Add 0.2 mL of MEM (+ 10% FBS)

↓

Cell count measurement

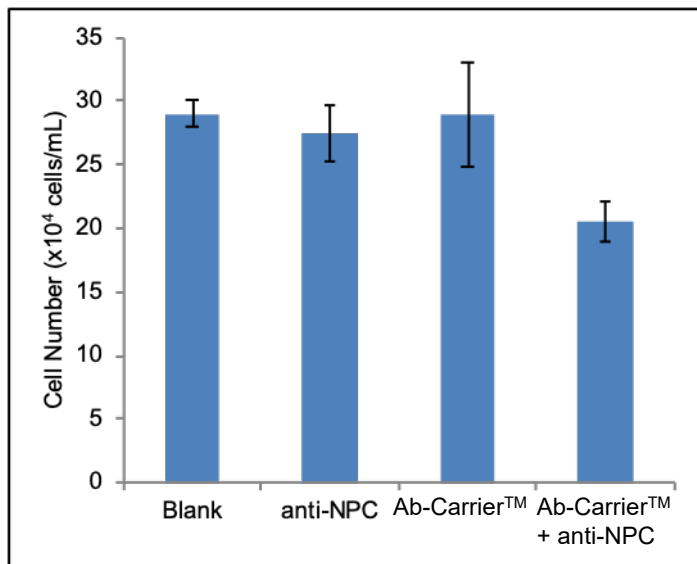


Figure 1. Inhibition of cell growth by anti-NPC antibody (n = 4)

Blank: unprocessed

anti-NPC: anti-NPC antibody only

Ab-Carrier™: only Ab-Carrier

Ab-Carrier™ + anti-NPC: mixture of Ab-Carrier™ and anti-NPC antibody

Antibodies used:

Anti-NPC antibody (Sigma-Aldrich)

Monoclonal Anti-Nuclear Pore Complex

Proteins, clone 414 (mouse IgG1)

The antibody was used after removing the preservative (NaN₃) by desalting.

24 hours after seeding HeLa cells (5×10^4 cells/mL) in a 96-well plate, a pre-mixture of antibody transfection reagent Ab-Carrier™ and anti-nuclear pore complex (NPC) antibody were added and cells were cultured for another 24 hours before comparing cell number with and without treatment. Compared to untreated cells, cells treated with a mixture of Ab-Carrier™ and anti-NPC antibody exhibited ~69% suppression of proliferation. We conclude that the antibody introduced into cells bound to the nuclear pore complex and impaired proliferation by inhibiting substance transport across the nuclear membrane.

Protenova Co., Ltd. 〒769-2604

Employee: 1488, Nishimura, Higashikagawa, Kagawa

WEB TEL 0879-49-0702 / FAX 0879-49-0703

WEB PAGE <http://protenova.com>