





Polymer Detection system Mouse on Mouse V-Histofine® MOUSESTAIN KIT

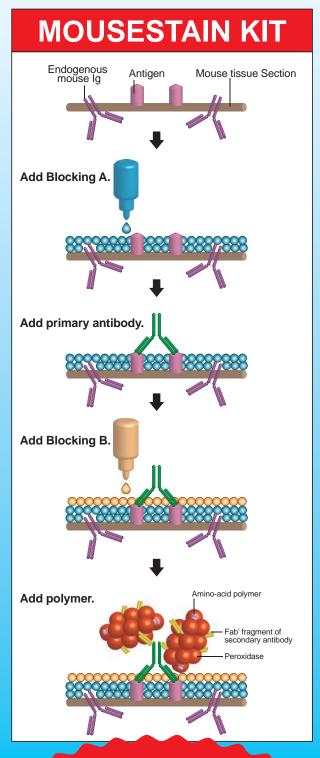

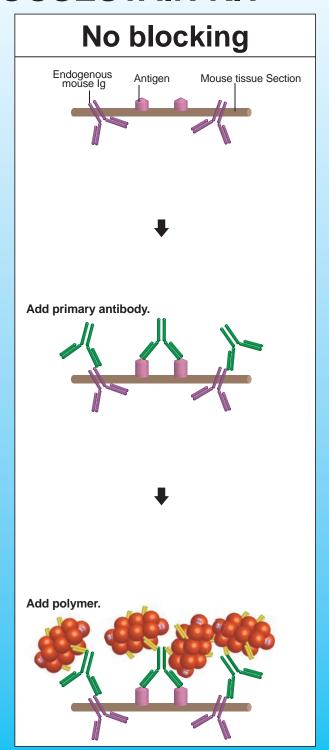

No background with double blocking


High background


Feature

✓-Histofine[®] MOUSESTAIN KIT is the mouse on mouse system. It is designed specifically to allow immunohistochemical staining with a mouse primary antibody on formalin-fixed paraffin-embedded mouse tissue sections. This kit consists of Blocking reagent A, Blocking reagent B, and Simple Stain Mouse Max PO (M) which is the labeled polymer prepared by combining amino acid polymer with multiple molecules of peroxidase and goat anti-mouse Ig which is reduced to Fab'. To eliminate background staining, this kit uses Blocking reagent A and Blocking reagent B.

Advantages


- 1. No reaction to endogenous mouse immunoglobulin in the tissue.
- 2. Simplified staining steps.
- 3. High sensitivity.
- 4. Unaffected by endogenous biotin in the tissue.
- 5. Ready to use.

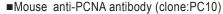

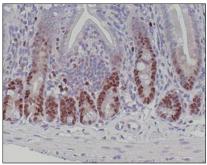

Mouse colon stained with **N**-Histofine® MOUSESTAIN KIT and DAB chromogen. Note nuclear staining of Epithelium cells.

■Mouse anti-Muscle Actin antibody (clone:HHF35)

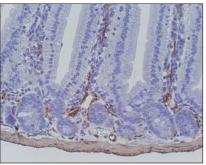

Mouse colon stained with **M**-Histofine® MOUSESTAIN KIT and DAB chromogen. Note intense staining of smooth muscle and muscularis propria in lamina propria mucosae.

Product List

Liquid. Ready to use.

Liquid. Ready to use					
	Product Name	Slides	Volume	Code	For use with
For Mouse Tissue Sections					
M -Histofine [®]	³ Simple Stain Mouse MAX PO (R)	170	1 x 17ml	414341F	Rabbit primary antibody
M -Histofine®	Simple Stain Mouse MAX PO (G)	170	1 x 17ml	414351F	Goat primary antibody
M -Histofine [®]	Simple Stain Mouse MAX PO (Rat)	170	1 x 17ml	414311F	Rat primary antibody
M _Histofine®	MOUSESTAIN KIT	50	3 x 6ml	414321F	Mouse
AW-I IISLUIIIIE		500	9 x 17ml	414322F	primary antibody
	Blocking Reagent A Blocking Reagent B Simple Stain Mouse Max PO (M)				
For Rat Tissue S	Sections				
N -Histofine®	Simple Stain Rat MAX PO (MULTI)	170	1 x 17ml	414191F	Mouse and rabbit primary antibodies
M -Histofine®	Simple Stain Rat MAX PO (M)	170	1 x 17ml	414171F	Mouse primary antibody
M -Histofine®	Simple Stain Rat MAX PO (R)	170	1 x 17ml	414181F	Rabbit primary antibody
M -Histofine [®]	Simple Stain Rat MAX PO (G)	170	1 x 17ml	414331F	Goat primary antibody
Enzyme Substra	ite Systems			,	
M -Histofine [®]	DAB-3S kit	500	3 x 3ml	415192F	Peroxidase
Reagent A : DAB solution concentrate Reagent B : Enhancer solution concentrate Reagent C : Hydrogen peroxide solution concentrate					
M -Histofine [®]	[®] Simple Stain AEC Solution		3 x 17ml 9 x 17ml	415182F 415184F	Peroxidase



Distributed in North America by:

Cosmo Bio USA

2792 Loker Ave W, Suite 101 Carlsbad, CA 92010

TEL: 760-431-4600 FAX: 760-431-4604 email: info@cosmobiousa.com web: www.cosmobiousa.com **NICHIREI BIOSCIENCES INC.**

6-19-20, TSUKIJI, CHUO-KU, TOKYO, 104-8402, JAPAN PHONE: 81-3-3248-2207 FAX: 81-3-3248-2243 http://www.nichirei.co.jp/bio/english

2016.07